HEXANIOS G+R IP sterile PAE

- Sans aldéhyde
- Produit biocide à large spectre
- Activités conformes aux normes européennes
- Testé sur surfaces et souches de terrain
- Conditionné en double emballage
- Radiostérilisé
- Aldehyde free
- Biocidal product with broad spectrum
- Activity in compliance with European standards
- Tested on field strains and critical surfaces
- Double-bagged packaged
- Gamma irradiated

INDICATIONS

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INDICATIONS

Cleaning and disinfection of floors, walls and surfaces in the Pharmaceutical Industry for clean rooms (A and B grades).

COMPOSITION

Polyhexamethylen biguanide hydrochloride (CAS N° 27083-27-8: 0.10 mg/g), didecyldimethylammonium chloride (CAS N°. 7173-51-5: 0.98 mg/g), excipients.

MICROBIOLOGICAL PROPERTIES

Bactericide in 5 min, 20°C (EN 1040, EN 1276, EN 13697).
Yeasticide in 15 min, 20°C (EN 1275, EN 1650, EN 13697).
Active at 20°C against BVDV (surrogate of HCV) and Herpes virus in 5 min and against HIV-1 and HBV in 10 min.

Biocidal activity on field surfaces and strains (bacteria, yeasts, moulds) according to EN 13697.

Other data concerning the product (material compatibilities...): see scientific file.

For each batch, Certificate of gamma irradiation, Certificate of analysis, LAL Testing analysis and Sterility report are available.

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PHYSICO-CHEMICAL DATA

- Clear solution
- Density at 20°C: approx 1
- Non-corrosive towards metals (non oxidising)
- pH of pure product: 7 to 8

INSTRUCTIONS FOR USE

Used undiluted.

Remove the first layer of packaging before entering into the clean room. Remove the second layer of packaging on first use.

Apply on the surface to be treated respecting the indicated contact time for the required antimicrobial activity. Apply in sufficient amount (30 to 40 ml/m²) and ensure that the surfaces remain wet throughout.

Rinse after use.

For frequency of use and cleaning of the application equipment, please refer to the hygiene plan in place.

PRECAUTIONS FOR USE

Dangerous. Follow the instructions for use (drawn up according to the European rules in force regarding the classification and labelling of chemical products). Use the biocides with precaution. Before use, read the label and the product information.

Store between +5°C and +35°C.

Biocidal product designed for surface disinfection (Group 1 - PT2) - Reserved for professional use.

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6 x 1L spray gamma irradiated double-bagged: 1139.534