SURFANIOS IP sterile PAE

- Biocide à large spectre PAE
- Association synergique de 2 principes actifs antimicrobiens
- Sans aldéhyde
- Activités conformes aux normes européennes
- Fabriqué en eau WFI
- Conditionné en double emballage
- Radiostérilisé
- Ready to use biocidal product with broad spectrum
- Synergistic combination of 2 antimicrobial active ingredients
- Aldehyde free
- Activity in compliance with European standards
- Made with WFI
- Double-bagged packaged
- Gamma irradiated

INDICATIONS

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INDICATIONS

Cleaning and disinfection of floors, walls, surfaces and equipment in the Pharmaceutical Industry for clean rooms (A and B grades).

COMPOSITION

N-(3-aminopropyl)-N-dodecylpropan-1,3-diamine - N° CAS 2372-82-9 (0,51 mg/g), Didecyldimethylammonium chloride (N° CAS 7173-51-5 (0,25 mg/g).

MICROBIOLOGICAL PROPERTIES

Bactericide in 5 min, 20°C (EN 1276, EN 13697).
Yeasticide in 5 min, 20°C (EN 1650).
Active in 5 min, 20°C against HIV-1, PRV (surrogate of HBV), BVDV (surrogate of HCV) and H1N1.

Biocidal activity on field surfaces and strains (bacteria, yeasts, moulds) according to EN 13697.

Other data concerning the product (material compatibilities...) : see scientific file.

For each batch, Certificate of gamma irradiation, Certificate of analysis, LAL Testing analysis and Sterility report are available.

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PHYSICO-CHEMICAL DATA

- Density at +20°C: approx. 1
- pH of product at 20°C: 9 to 12
- Non corrosive (free of oxidising agents)

INSTRUCTIONS FOR USE

To be used undiluted.

Remove the first package before entering into the clean room. Remove the second package on first use. Apply in sufficient amount (+/- 30 ml/m²) respecting the indicated contact time, for the required antimicrobial activity. If necessary remove the surplus after the contact time. For frequency of use and cleaning of the application equipment, please refer to the hygiene plan in place.

PRECAUTIONS FOR USE

Dangerous - Please follow the instructions for use (drawn up according to the European rules in force regarding the classification and labelling of chemical products). Use the biocides with precaution. Before use, read the label and the product information.
Store between +5°C and +35°C.
Biocidal product designed for surface disinfection (Group 1 - PT2) - Reserved for professional use.

CONDITIONNEMENTS/PACKAGING

- 4 x 5L double emballage radiostérilisé : 1342.278
- 4 x 5L gamma irradiated double-bagged: 1342.278
- 6 x 1L double emballage radiostérilisé : 1342.534
- 6 x 1L gamma irradiated double-bagged: 1342.534