BACTERANIOS SF IP sterile concentré

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- *Biocidal product with broad spectrum*
- *Synergistic combination of 2 antimicrobial active ingredients*
- *Formaldehyde free*
- *Activity in compliance with European standards*
- *Double-bagged packaged*
- *Gamma irradiated*

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INDICATIONS

Cleaning and disinfection of floors, walls, surfaces and equipment in the Pharmaceutical Industry for clean rooms (A and B grades).

COMPOSITION

Glutaraldehyde (CAS N°. 111-30-8 : 50 mg/g), didecyldimethylammonium chloride (CAS N°. 7173-51-5: 120 mg/g).

MICROBIOLOGICAL PROPERTIES

Bactericide in 5 min, 0.5%, 20°C (EN 1276, EN 13697).
Yeasticide in 15 min, 0.5%, 20°C (EN 1650, EN 13697).
Fungicide in 30 min, 2%, 20°C (EN 1650, EN 13697).
Active in 5 min 0.5% 20°C against BVDV (surrogate of HCV) and PRV (surrogate of HBV).

Biocidal activity on field surfaces and strains (bacteria, yeasts, moulds) according to EN 13697.

Other data concerning the product (material compatibilities...): see scientific file.

For each batch, Certificate of gamma irradiation, Certificate of analysis, LAL Testing analysis and Sterility report are available.

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PHYSICO-CHEMICAL DATA

- Liquid colourless solution
- Soluble in all proportions in water
- Density at 20°C : approx. 1
- pH of product at 20°C : approx. 4
- Non-corrosive towards metals (non oxidising).

INSTRUCTIONS FOR USE

Concentrated solution.

Remove the first package before entering into the clean room. Remove the second package on first use.

Use at dilution from 0,5% (5ml/l) to 2% (20ml/l), i.e. 1 to 4 bottles for 12l of cold water, according to the required antimicrobial activity, respecting the indicated contact time. Apply a sufficient amount (+/- 30 ml/m²).

If necessary, remove the excess after the contact time. For frequency of use and cleaning of the application equipment, please refer to the hygiene plan in place.

PRECAUTIONS FOR USE

Dangerous. Follow the instructions for use (in accordance with Directive 99/45/EC and its adapted versions). Use the biocides with precaution. Before use, read the label and the product information.

Store between +5°C and +35°C.

Biocidal product designed for surface disinfection (Group 1 - PT2) - Reserved for professional use.

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