SURFANIOS PREMIUM IP sterile PAE

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INDICATIONS

Cleaning and disinfection of floors, walls and equipment in the Pharmaceutical Industry for clean rooms (A and B grades).

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Bactericide in 5 min, 20°C (EN 1276, EN 13697).

Yeasticide in 15 min, 20°C (EN 1650, EN 13697).

Active against PRV (surrogate of HBV), BVDV (surrogate of HCV), HIV-1, Herpesvirus, Vaccinia virus, Rotavirus in 5 min, 20°C; Coronavirus in 15 min, 20°C.

Other data concerning the product (material compatibilities...) : see scientific file.

For each batch, Certificate of gamma irradiation, Certificate of analysis, LAL Testing analysis and Sterility report are available.

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PHYSICO-CHEMICAL DATA

- pH: approx. 10.5
- Density: approx. 1.00
- Refraction index: approx. 1.33

INSTRUCTIONS FOR USE

To be used undiluted.

Remove the first package before entering into the clean room.

Remove the second package on first use.

Apply on the surface to be treated.

Respect the indicated contact time for the required antimicrobial activity.

Apply in sufficient amount (+/- 30 ml/m²).

If necessary remove the surplus after the contact time.

For frequency of use and cleaning of the application equipment, please refer to the hygiene plan in place.

PRECAUTIONS FOR USE

Dangerous - follow the instructions for use (drawn up according to the European rules in force regarding the classification and labelling of chemical products). Use the biocides with precaution. Before use, read the label and the product information.

Store between +5°C and +35°C.

Biocidal product designed for surface disinfection (Group 1 - PT 2) - Reserved for professional use.

CONDITIONNEMENTS/PACKAGING

6 flacons de 1L, double emballage radiostérilisé : 2409.534
6 bottles, 1L each, gamma irradiated double-bagged: 2409.534

6 flacons de 1L adaptable sur balai, double emballage radiostérilisé : 2409.559
6 bottles broom suitable, 1L each, gamma irradiated double-bagged: 2409.559

4 x 5L double emballage radiostérilisé : 2409.278
4 x 5L gamma irradiated double-bagged: 2409.278