SURFANIOS PREMIUM IP STERILE PAE

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INDICATIONS

Cleaning and disinfection of floors, walls and equipment in the Pharmaceutical Industry for clean rooms (A and B grades).

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Active against PRV (surrogate of HBV), BVDV (surrogate of HCV), HIV-1, Herpesvirus, Vaccinia virus, Rotavirus in 5 min, 20°C; Coronavirus in 15 min, 20°C.

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Other data concerning the product (material compatibilities...) : see scientific file.

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For each batch, Certificate of gamma irradiation, Certificate of analysis, LAL Testing analysis and Sterility report are available.

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PHYSICO-CHEMICAL DATA

- pH: approx. 10.5
- Density: approx. 1.00
- Refraction index: approx. 1.33

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INSTRUCTIONS FOR USE

To be used undiluted.

Remove the first package before entering into the clean room.

Remove the second package on first use.

Apply on the surface to be treated.

Respect the indicated contact time for the required anti-microbial activity.

Apply in sufficient amount (+/- 30 ml/m²).

If necessary remove the surplus after the contact time.

For frequency of use and cleaning of the application equipment, please refer to the hygiene plan in place.

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PRECAUTIONS FOR USE

Dangerous - follow the instructions for use (drawn up according to the European rules in force regarding the classification and labelling of chemical products). Use the biocides with precaution. Before use, read the label and the product information.

Store between +5°C and +35°C.

Biocidal product designed for surface disinfection (Group 1 - PT 2) - Reserved for professional use.

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6 bottles, 1L each, gamma irradiated double-bagged: 2409.534

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6 bottles broom suitable, 1L each, gamma irradiated double-bagged: 2409.559

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4 x 5L gamma irradiated double-bagged: 2409.278

Summary

Product Brochure	Page	7
Regulation and General Information Brochure	Page	9
Composition	Page	11
Stability and storage conditions	Page	13
Studies and expertises		
- Microbiology	Page	15
- Compatibilities	Page	21
Regulation	Page	23

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***SURFANIOS PREMIUM IP sterile PAE** is the ready to use form of SURFANIOS PREMIUM IP sterile.*

*Consequently, microbiological properties of SURFANIOS PREMIUM IP sterile are applicable to **SURFANIOS PREMIUM IP sterile PAE** and reciprocally.*

***SURFANIOS PREMIUM IP sterile PAE** is filtered and packaged in clean rooms, and terminally gamma irradiated.*

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Regulation and general information

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***SURFANIOS PREMIUM IP STERILE PAE** is designed, produced and controlled by the Laboratoires ANIOS, certified by the AFAQ Organism under the number 1995/3723 in accordance with the ISO 9001 Quality System*

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***SURFANIOS PREMIUM IP STERILE PAE** is in compliance with the requirements of the European regulation related to biocides for disinfection of surfaces (Group 1 – PT 2)*

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***SURFANIOS PREMIUM IP STERILE PAE** formulation fulfils the requirements of Regulation (EC) No 648/2004 on detergents and its amendments*

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***SURFANIOS PREMIUM IP STERILE PAE** is labelled in accordance with the European regulation related to the classification and labelling of chemical products*

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***SURFANIOS PREMIUM IP STERILE PAE** meets our voluntary eco-design commitment included in our ANIOSAFE charter*

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SURFANIOS PREMIUM IP

Stérile PAE

Stabilité et conditions de conservation (1)
Stability and storage conditions (1)

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SURFANIOS PREMIUM IP

Stérile PAE

Stabilité et conditions de conservation (2) *Stability and storage conditions (2)*

Pure product ready to use :

- Storage between +5°C and +35°C.

- Physico-chemical stability:

The shelf-life of SURFANIOS PREMIUM IP Sterile PAE was determined on the basis of physico-chemical stability studies and active substances titration in recommended storage conditions.

- Not opened can of 1L and 5L: 2 years from the manufacturing date indicated on the label.

- Opened can of 1L and 5L : 3 months after opening

- Microbiological stability:

- Opened can of 1L and 5L: the sterility of the contents after opening must be evaluated by the user according to its terms of use.

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*SURFANIOS PREMIUM IP STERILE PAE is the ready to use form of
SURFANIOS PREMIUM IP STERILE CONCENTRE at the concentration of $1,2 \pm 0,2\%$*

Etudes sur produit concentré : <i>Studies on concentrated product :</i> SURFANIOS PREMIUM – SURFANIOS PREMIUM IP stérile	Résultats <i>Results</i>		
	Concentration active <i>Active concentration</i>	Temps de contact <i>Contact time</i>	Conditions spécifiques <i>Specific conditions</i>
BACTERICIDIE <i>Bactericidal activity</i>			
EN 1040	0,10 %	5 min.	
EN 1276	0,25 %	5 min.	Conditions de saleté <i>Dirty conditions</i>
EN 13697	1,00 %	5 min.	Conditions de saleté
	0,50 %	15 min.	<i>Dirty conditions</i>
	0,25 %	30 min.	
<u>Souches additionnelles</u> <u>Additional strains</u>			
Acinetobacter baumannii BLSE selon / according to EN 1276	0,10 %	5 min.	Conditions de saleté <i>Dirty conditions</i>
Enterobacter aerogenes BLSE selon / according to EN 13697	0,10 %	15 min.	Conditions de saleté <i>Dirty conditions</i>
Enterobacter cloacae BLSE selon / according to EN 13697	0,25 %	15 min.	Conditions de saleté <i>Dirty conditions</i>
Enterococcus faecium ERV / VRE selon / according to EN 1276	0,10 %	5 min.	Conditions de saleté <i>Dirty conditions</i>
Escherichia coli BLSE selon / according to EN 13697	0,25 %	15 min.	Conditions de saleté <i>Dirty conditions</i>
Klebsiella pneumoniae BLSE selon / according to EN 13697	0,25 %	15 min.	Conditions de saleté <i>Dirty conditions</i>
Staphylococcus aureus SARM / MRSA selon / according to EN 1276	0,05 %	5 min.	Conditions de saleté <i>Dirty conditions</i>

Efficacité antimicrobienne – SURFANIOS PREMIUM IP stérile PAE - Page 1/3 - le 11.04.2016
Antimicrobial effectiveness - SURFANIOS PREMIUM IP stérile PAE - Page 1/3 – on 11.04.2016

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*SURFANIOS PREMIUM IP STERILE PAE is the ready to use form of
 SURFANIOS PREMIUM IP STERILE CONCENTRE at the concentration of $1,2 \pm 0,2\%$*

Etudes sur produit concentré : <i>Studies on concentrated product :</i> SURFANIOS PREMIUM – SURFANIOS PREMIUM IP stérile	Résultats <i>Results</i>		
	Concentration active <i>Active concentration</i>	Temps de contact <i>Contact time</i>	Conditions spécifiques <i>Specific conditions</i>
TUBERCULOCIDIE <i>Tuberculocidal activity</i>			
EN 14348 : Mycobacterium terrae	0,25 %	60 min.	Conditions de saleté
	0,50 %	30 min.	<i>Dirty conditions</i>
EN 14563 : Mycobacterium terrae	0,25 %	60 min.	Conditions de saleté
	0,50 %	30 min.	<i>Dirty conditions</i>
LEVURICIDIE et/ou FONGICIDIE <i>Yeasticidal and/or fungicidal activity</i>			
EN 1275 : Candida albicans Aspergillus niger	0,05 % 0,25 %	15 min. 15 min.	Conditions de saleté <i>Dirty conditions</i> Conditions de saleté <i>Dirty conditions</i>
EN 1650 : Candida albicans Aspergillus niger	0,25 % 0,25 %	15 min. 120 min.	
EN 13697 : Candida albicans	0,50 % 0,25 %	15 min. 30 min.	
Aspergillus niger	1,00 % 0,25 %	15 min. 30 min.	

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Not concerned by biodegradability notion

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0,2 %¹

*Readily biodegradable organic material*²

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Biodegradability within 28 days or less:

- ≥ 60% according to respirometric methods³ (assessment of CO₂ production or oxygen consumption)
- ≥ 70% according to methods measuring dissolved organic carbon³ (DOC) (assessment of DOC die away)

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Substances which not fulfil the readily biodegradability criteria (but possibly displaying inherent biodegradability criteria⁴), or substances without any biodegradability data.

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SURFANIOS PREMIUM IP sterile PAE contains 100 % of inorganic material and of readily biodegradable organic material.

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HAZARDS IDENTIFICATION

Health

*This mixture does not present a health hazard with the exception of possible occupational exposure thresholds.**

Environment

*This mixture does not present an environmental hazard. No known or foreseeable environmental damage under standard conditions of use.**

* Classification drawn up in accordance with European regulation related to the classification and labelling of chemical products.

PERSONAL PROTECTION UNDER THE CONDITIONS OF THE INTENDED USE

Refers exclusively to this product. If the product is handled and/or exposed simultaneously with other chemical agents, these must be taken into consideration when choosing personal safety equipment.

Respiratory protection

Not applicable in normal conditions of use.

Hand protection

When handling this product, wear suitable gloves (nitrile, latex or vinyle) impervious in accordance with standard EN374

Gloves must be replaced immediately if they show signs of wear and tear

Eye and face protection

Avoid contact with eyes

Water source nearby

*** please refer to SDS §8 for TWA (Time-Weighted Average)/STEL (Short-term exposure limit) of concerned substances present in the product*

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The Safety Data Sheet (SDS) of this product is available on our Web site www.anios.com. In addition to visualization and/or impression, it is possible for you to download it.

The consultation of this SDS via internet enables you to receive later on by email any updated version of this document.

In the absence of an Internet access, the SDS remains available on request near our laboratories (tel.: +33 3 20 67 67 67 - fax: +33 3 20 67 67 68)

* an access is also possible at the site <http://www.quickfds.com>