SURFANIOS IP sterile concentré

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- *Aldehyde free*
- *Biocidal product with broad spectrum*
- *Activity in compliance with European standards*
- *Tested on field strains and critical surfaces*
- *Double-bagged packaged*
- *Gamma irradiated*

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CHARACTERISTICS

Cleaning and disinfection of floors, surfaces, walls and equipment in the Pharmaceutical Industry (Clean rooms).

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Delivered, for each batch, with Certificate of gamma irradiation, Certificate of analysis, LAL Testing analysis and Sterility report.

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PHYSICO-CHEMICAL DATA

- Density at +20°C : approx. 1
- pH of pur product at 20°C : > 11
- pH of product at 0.25% : approx. 8.5 (in hard water)
- Non corrosive (non oxidising)

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INSTRUCTIONS FOR USE

Concentrated solution.

Remove the first layer of packaging before entering into the clean room. Remove the second layer of packaging on first use.

Use at dilution of 0,25% (2,5ml/l) to 1% (10ml/l) according to the required antimicrobial activity, respecting the indicated contact time.

Apply in sufficient amount (30 to 40 ml/m²) and ensure that the surfaces remain wet throughout.

Rinse after use.

For frequency of use and cleaning of the application equipment, please refer to the hygiene plan in place.

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PRECAUTIONS FOR USE

Dangerous - follow the instructions for use (in accordance with Directive 99/45/EC and its adapted versions). Use the biocides with precaution. Before use, read the label and the product information.

Store between +5°C and +35°C.

Biocidal product designed for surface disinfection (Group 1 - PT2 - Directive 98/8/EC) - Reserved for professional use

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15 x 8 doses of 15 ml gamma irradiated double-bagged: 871.245