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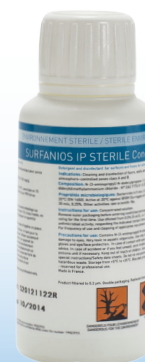
INDICATIONS

Cleaning and disinfection of floors, surfaces, walls and equipment in the Pharmaceutical Industry for clean rooms (A and B grades).

COMPOSITION

N-(3-aminopropyl)-N-dodecylpropan-1,3-diamine – N° CAS 2372-82-9 (51 mg/g), Didecyldimethylammonium chloride (N° CAS 7173-51-5 (25 mg/g).

MICROBIOLOGICAL PROPERTIES

Bactericide in 5 min, 0,5% 20°C (EN 1276, EN 13697).
Yeasticide in 5 min, 0,25%, 20°C (EN 1650).
Active in 5 min, 0,25%, 20°C against HIV-1, PRV (surrogate of HBV), BVDV (surrogate of HCV) and H1N1.

Biocidal activity on field surfaces and strains (bacteria, yeasts, moulds) according to EN 13697.

Other data concerning the product (material compatibilities...) : see scientific file.

For each batch, Certificate of gamma irradiation, Certificate of analysis, LAL Testing analysis and Sterility report are available.

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PHYSICO-CHEMICAL DATA

- Density at +20°C: approx. 1
- pH of pure product at 20°C: > 11
- Non corrosive (non oxidising)

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INSTRUCTIONS FOR USE

Concentrated solution.

Remove the first package before entering into the clean room. Remove the second package on first use. Use at dilution of 0,5% (5ml/l), i.e. 1 flask for 12L of cold water, respecting the indicated contact time, for the required antimicrobial activity. Apply in sufficient amount (+/- 30 ml/m²). If necessary remove the surplus after the contact time. For frequency of use and cleaning of the application equipment, please refer to the hygiene plan in place.

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PRECAUTIONS FOR USE

Dangerous - Please follow the instructions for use (drawn up according to the European rules in force regarding the classification and labelling of chemical products).

Use the biocides with precaution. Before use, read the label and the product information.

Store between +5°C and +35°C.

Biocidal product designed for surface disinfection (Group 1 - PT2) - Reserved for professional use.

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