

Premier Klerwater-WFI Sterile WFI Quality Water

Bulk quantities of sterile spore free WFI quality water for residue removal and dilution of sterile biocide concentrates.

Advantages

■ Suitable for use in product contact areas

Premier Klerwater-WFI Sterile WFI Quality Water is produced from the highest grade (EP – which requires a distillation process) Water for Injection, with guaranteed endotoxin levels below 0.25 EU/ml so is suitable for use in product contact areas. It is double bagged for easy transfer into Grade A and B cleanrooms.

■ Suitable for dilution of sterile biocide concentrates

Klerwater-WFI is useful for the dilution of sterile concentrates to maintain the required quality of diluted product for high grade cleanrooms.

■ Useful for rinsing to remove residues

Klerwater-WFI is an effective way of eliminating any residues after disinfection. Wipe down with Klerwater-WFI after using the Klercide range of biocides in areas where residues may be an issue. Klerwater-WFI acts as an excellent solvent, it is particularly effective at removing ionic and polar covalent compounds.

■ Sterile and quality assured

The product is manufactured to a tight specification. Gamma irradiation is carried out to a validated procedure. Samples from each batch are tested for sterility and a quantitative endotoxin analysis is carried out before QC release.

■ Cost effective

The sterile water is available in either 1 litre trigger sprays fitted with the SteriShield Delivery System (SDS) and bulk 5 litre capped bottles. The SDS ensures all the contents are dispensed from the bottle so there is no wastage. The SDS has been validated to give an in-use shelf life of three months so there is no need to discard unused product.



Description

Premier Klerwater-WFI Sterile WFI Quality Water is 0.2 micron filtered, filled and double bagged in a Grade C (ISO Class 7) cleanroom before irradiation using a validated process at no less than 25 kGy to ensure it is spore free.

The 1 litre trigger sprays incorporate the patented Sterishield Delivery System and an adjustable trigger spray. The unique way in which the SDS works as a closed system ensures that the sterility of the contents is preserved throughout use. Also available in 5 litre capped bottles for dilution of sterile biocide concentrates.

Validation

Manufactured by an ISO 9001:2008, ISO 13485:2003 and European Medical Device Directive 93/42/EEC accredited company.

Certificates of analysis, including quantitative endotoxin levels <0.25 EU/ml, irradiation available with every batch. Sterility and endotoxin tested to an approved pharmacopoeial method.

2 year shelf life.

Product Codes

CODE	OLD CODE	DESCRIPTION	SIZE	UNIT OF SALE
3056490		Premier Klerwater-WFI Sterile WFI Quality Water Spray	1.0 litre	6 bottles x 1.0L
3056500		Premier Klerwater-WFI Sterile WFI Quality Water Capped	5.0 litre	4 bottles x 5.0L

For more information please contact info@shieldmedicare.com or visit www.shieldmedicare.com