



T +44 (0) 2920 854 390  
F +44 (0) 2920 854 391

BRUNEL WAY, BAGLAN ENERGY PARK,  
NEATH, SA11 2GA, UK  
www.ecolabcc.com  
info@ecolab.com

# Klercide™

## 70/30 Sterile Alcohol

**Sterile spore free 70/30 v/v alcohol, either IPA or Denatured Ethanol blended with deionised water.**

### Advantages

#### ■ Easy to use

Convenient range of comfortable and easy to use trigger sprays is available, with lightweight options from 500 ml to 4 litres. The large droplet size reduces the risk of inhalation.

#### ■ Cost effective

All the sterile alcohol can be dispensed from the bottle, so there is no wastage. The SteriShield Delivery System has been validated to give an in-use shelf life of three months, so there is no need to discard unused product.

#### ■ Optimum surface coverage

The fully adjustable trigger spray allows the liquid to be dispensed as either a jet or spray, enabling thorough wetting of the surface so that disinfection procedures are effective.

#### ■ More environmentally friendly

The product contains no propellants so does not require any special disposal procedures. Using a 4 litre spray significantly reduces the packaging waste.



### Description

Klercide 70/30 IPA is a blend of 70% v/v Isopropyl Alcohol with deionised water. Klercide 70/30 Denatured Ethanol is 70% v/v denatured ethanol blended with deionised water.

The alcohol blend is 0.2 micron filtered, filled and double bagged in a Grade C (ISO class 7) cleanroom, before irradiation using a validated process at no less than 25 kGy to ensure a spore free alcohol.

All trigger spray bottles incorporate the patented SteriShield Delivery System and an adjustable trigger spray. The unique way in which the SteriShield Delivery System works as a closed system, ensures that the sterility of the contents is preserved throughout use.

## Protocol

Ready to use, so can be sprayed directly onto the surface to be sanitised. Always ensure the whole surface is in contact with the liquid. For optimum results, wipe the surface to physically remove any contaminants.

*Use biocides safely. Always read the label and product information before use.*

## Validation

Manufactured by a BS EN ISO 9001:2008, BS EN ISO 13485:2012 and European Medical Device Directive 93/42/EEC accredited company.

Passes CEN test 1276:1997 for bactericidal efficacy.

Certificates of analysis, irradiation and sterility with every batch. Sterility tested to a method approved by the United States Pharmacopœia.

2 year shelf life.

In use sterility guaranteed by the SteriShield Delivery System, validated for three months.

Full supporting documentation is available on these products.

## Product Codes

| CODE    | OLD CODE | LANGUAGES ON LABEL | DESCRIPTION                      | WATER QUALITY | STERILE / FILTERED | SIZE      | UNIT OF SALE      |
|---------|----------|--------------------|----------------------------------|---------------|--------------------|-----------|-------------------|
| 3078260 | 3035330  | GB;DE;FR;IT;ES;NL  | Klercide 70/30 Denatured Ethanol | DI            | Sterile            | 500ml     | 12 bottles x 0.5L |
| 3078270 | 3055600  | GB;SE;NO;FI;PL;DK  | Klercide 70/30 Denatured Ethanol | DI            | Sterile            | 500ml     | 12 bottles x 0.5L |
| 3078240 | 3055610  | GB;SE;NO;FI;PL;DK  | Klercide 70/30 Denatured Ethanol | DI            | Sterile            | 1.0 litre | 6 bottles x 1.0L  |
| 3078280 | 3035360  | GB;DE;FR;IT;ES;NL  | Klercide 70/30 Denatured Ethanol | DI            | Sterile            | 5.0 litre | 4 bottles x 5.0L  |
| 3078290 | 3035350  | GB;DE;FR;IT;ES;NL  | Klercide 70/30 Denatured Ethanol | DI            | Sterile            | 4.0 litre | 4 bottles x 4.0L  |
| 3078300 | 3035340  | GB;DE;FR;IT;ES;NL  | Klercide 70/30 Denatured Ethanol | DI            | Sterile            | 1.0 litre | 6 bottles x 1.0L  |
| CODE    | OLD CODE | LANGUAGES ON LABEL | DESCRIPTION                      | WATER QUALITY | STERILE / FILTERED | SIZE      | UNIT OF SALE      |
| 3078440 | 3035290  | GB;DE;FR;IT;ES;NL  | Klercide 70/30 IPA               | DI            | Sterile            | 500ml     | 12 bottles x 0.5L |
| 3078450 | 3055560  | GB;SE;NO;FI;PL;DK  | Klercide 70/30 IPA               | DI            | Sterile            | 1.0 litre | 6 bottles x 1.0L  |
| 3078460 | 3035320  | GB;DE;FR;IT;ES;NL  | Klercide 70/30 IPA               | DI            | Sterile            | 5.0 litre | 4 bottles x 5.0L  |
| 3078470 | 3035310  | GB;DE;FR;IT;ES;NL  | Klercide 70/30 IPA               | DI            | Sterile            | 4.0 litre | 4 bottles x 4.0L  |
| 3078480 | 3035300  | GB;DE;FR;IT;ES;NL  | Klercide 70/30 IPA               | DI            | Sterile            | 1.0 litre | 6 bottles x 1.0L  |

*For more information please contact [info@ecolab.com](mailto:info@ecolab.com) or visit [www.ecolabcc.com](http://www.ecolabcc.com)*