

Versatile, High Grade Disinfection with Easy Application



Klercide™ 70|30 Alcohol Range Blended with WFI

Sterile, spore free 70|30 Alcohol blended with Water for Injection (WFI)

A choice of three blends of IPA, Denatured Ethanol or Pharma Ethanol, with the highest grade (EP) Water for Injection providing a versatile, high-quality product with guaranteed endotoxin levels below 0.25 EU/ml, for use in Grade A and B cleanrooms.

Available in a range of formats for optimum performance depending on the application method required.

The Klercide 70|30 Alcohol blended with WFI range is:

- ▲ Triple bagged for easy transfer into Grade A and B cleanrooms
- ▲ Gamma irradiated to ensure a sterile, spore free product
- ▲ Endotoxin tested to an approved pharmacopoeial method
- ▲ Fitted with the SteriShield Delivery System on the trigger spray to give an in-use shelf life of three months
- ▲ Available in a range of formats





Klercide™ 70|30 Alcohol Range Blended with WFI

Sterile, spore free 70|30 Alcohol blended with WFI

Why use the Klercide 70|30 Alcohol Range?

EFFECTIVE

Blended with WFI, Klercide 70|30 IPA is 70% isopropyl alcohol, Klercide 70|30 Denatured Ethanol is 70% denatured ethanol and Klercide 70|30 Pharma Ethanol is 70% ethanol denatured with IPA for a very low residue product.

OPTIMUM SURFACE COVERAGE

The choice of formats allows the user to select the most appropriate method of application, from the fine mist of an aerosol*, to a fully adjustable trigger spray which allows the liquid to be dispensed as either a jet or spray.

TOTAL RELIABILITY

The alcohol blend is 0.2 micron filtered and filled under Grade A laminar flow and bagged in a Grade C cleanroom before irradiation using a validated process at no less than 25 kGy. The trigger spray has the option of a large droplet size to reduce the risk of inhalation.

ECONOMICAL

All trigger spray bottles incorporate the patented SteriShield Delivery System (SDS) and an adjustable trigger spray. The SDS's unique closed system, ensures that the sterility of the contents is preserved throughout use and there is no wastage.



Klercide 70|30 IPA Range Blended with WFI
360ml Aerosol, 500ml, SDS Spray
1L SDS Spray, 5L Capped

VALIDATION

Manufactured by an ISO 9001:2008 and ISO 14001:2015 accredited company, the Klercide Alcohol range provides bactericidal, fungicidal and yeasticidal efficacy.

Certificates of Analysis and Conformity with quantitative endotoxin levels <0.25 EU/ml are supplied with every batch and the sterility test method is to an approved pharmacopoeial method.

With a two year shelf life (for triggers and capped product) and three years for aerosols, the products' in use sterility, is guaranteed by the SteriShield Delivery System, which is validated to enable a three month in-use shelf life to be assigned.

Full supporting documentation is available on these products.

*Not available in Pharma Ethanol range



Klercide 70|30 Pharma Ethanol Range Blended with WFI
1L SDS Spray, 5L Capped



Klercide 70|30 Denatured Ethanol Range Blended with WFI
360ml Aerosol, 500ml SDS Spray
1L SDS Spray, 5L Capped

WORLDWIDE HEADQUARTERS

1 Ecolab Place
St. Paul, MN 55102
USA

www.ecolab.com/lifesciences

© 2018 Ecolab USA Inc. All rights reserved. 05JAN18/EU

EUROPE HEADQUARTERS

Richtistr. 7
8304 Wallisellen
Switzerland

www.ecolablifesciences.com

EFFICIENT.
POWERFUL.
ECONOMICAL.

APPLICATION AREA

For use on floors, walls and hard surfaces.

INSTRUCTIONS FOR USE

Aerosol: Hold the aerosol upright and spray from a distance of 10cm to 20cm.

Spray: For optimum results hold bottle upright and spray from a distance of 10cm to 20cm. Always close the nozzle after use.

Wetting: For optimum results ensure complete wetting of the surfaces using the spray and wipe technique (max. 40ml/m²).

INGREDIENTS

See individual products for details.

PACKAGING INFORMATION

UNIT SIZE	UNIT OF SALE
Aerosol 360ml	24 cans x 360ml
500ml Spray	12 bottles x 500ml
1.0 litre spray	6 bottles x 1.0L
5.0 litre capped	4 bottles x 5.0L

FOR FULL PRODUCT ORDER CODES, PLEASE SEE OUR PRODUCT CATALOGUE OR VISIT OUR WEBSITE

ALWAYS READ THE LABEL AND PRODUCT INFORMATION BEFORE USE.

