

# Rapid Disinfection of Surfaces, Equipment and Production Components



## Klerwipe™ Impregnated Sterile Biocide Wipes

Ergonomic, simple to use pouch wipes with a wide range of applications

These sterile, low-particulate pouch wipes are impregnated with either Klercide Quat|Biguanide, Sporicidal Chlorine|Quat, Sporicidal Low Residue Peroxide or Amine in a convenient resealable pouch.

They are highly absorbent, allowing rapid disinfection of surfaces, and they remain wet throughout use ensuring a consistent application of biocide to the whole surface area.

Klerwipe Biocide Wipes provide:

- ▲ Very effective biocides including one with sporicidal activity
- ▲ Effective rotational use
- ▲ Klertear easy open feature - ideal for gloved hands
- ▲ Easy to use, resealable pouches for longer life
- ▲ Ideal for sterile environments





# Klerwipe™ Impregnated Sterile Biocide Wipes

Ergonomic, simple to use pouch wipes with a wide range of applications

## Why use Klerwipe Impregnated Sterile Biocide Wipes?

### EFFECTIVE

All wipes are impregnated with 0.2 micron filtered fluids from the Klercide range of biocides.

They are double bagged for assured transfer and provide optimal effectiveness against a wide range of organisms commonly encountered in a cleanroom.

### LOW RISK

The selected biocides are non-toxic and non-corrosive under normal use, so can be used in most cleanroom applications.

Being pre-impregnated, Klerwipe pouch wipes eliminate the potential risk, when mixing a concentrate, of a dilution error, and deliver the correct concentration of active ingredient to the surface being disinfected.

The Klerwipe pouch wipes are either packed in a Grade C (ISO Class 7) cleanroom and sterilised by gamma irradiation or aseptically processed, under a Grade A (ISO Class 5) Laminar Air Flow system to a validated procedure.

### SIMPLE TO USE

The 200mm x 200mm wipes are made from a highly absorbent, non-woven, binder free, hydroentangled cellulose/ polyester blend.

Each pouch is designed for sessional use and contains 15 wipes so there is no wastage at the end of a manufacturing session. The pouch is easily resealable to prevent the wipes drying out. The pouch itself does not generate any particulate contamination.

### VALIDATION

Manufactured by a ISO 9001:2008 and ISO 14001:2015 accredited company.

Certificates of Analysis and Sterility are available for every batch of Klerwipe pouch wipes.

Klerwipe Quat / Biguanide and Amine - two year shelf life.

Klerwipe Sporicidal Chlorine / Quat and Sporicidal Low Residue Peroxide - one year shelf life.



Klerwipe Sterile Impregnated Biocide Pouch Wipes

### WORLDWIDE HEADQUARTERS

1 Ecolab Place  
St. Paul, MN 55102  
USA  
www.ecolab.com

### EUROPE HEADQUARTERS

Richtistr. 7  
8304 Wallisellen  
Switzerland  
www.ecolablifesciences.com

EFFECTIVE.  
SIMPLE.  
CONVENIENT.

### APPLICATION AREA

For use on hard surfaces.

### INSTRUCTIONS FOR USE

For optimum results ensure complete wetting of the surfaces. Re-seal when not in use. Dispose of packaging only if emptied and closed.

### INGREDIENTS

See individual products for details.

### PACKAGING INFORMATION

UNIT SIZE	UNIT OF SALE
200 x 200mm Wipe	20 packs x 15 wipes

FOR FULL PRODUCT ORDER CODES, PLEASE SEE OUR PRODUCT CATALOGUE OR VISIT OUR WEBSITE

USE BIOCIDES SAFELY. ALWAYS READ THE LABEL AND PRODUCT INFORMATION BEFORE USE.

