



T +44 (0) 2920 854 390
F +44 (0) 2920 854 391

BRUNEL WAY, BAGLAN ENERGY PARK,
NEATH, SA11 2GA, UK
www.ecolabcc.com
infocc@ecolab.com

Klercide™

Quat / Biguanide

A sterile broad spectrum rotational disinfectant effective even under conditions of heavy soiling.

Advantages

■ Very effective rotational biocide

Klercide Quat / Biguanide has a good broad spectrum of activity, which will control most microbial contamination likely to be found in a cleanroom environment. It can be used safely and effectively as part of a rotation with any other Ecolab Contamination Control biocide.

■ Easy to use

Klercide Quat / Biguanide is classified as non-irritant and non-corrosive so it can be used in most cleanroom applications. Ready to use products and concentrates are available for use in large facilities.

■ Cost effective

All the biocide can be dispensed from the bottle so there is no wastage.

■ Environmentally friendly

The packaging does not contain any propellant so does not require any special disposal procedures.



Description

Klercide Quat / Biguanide is a sterile cleanroom biocide. The active ingredients are a mix of a quaternary ammonium compound and a biguanide. The blend gives broad spectrum activity even under conditions of heavy organic soiling. No additional activators are required. Klercide Quat / Biguanide is suitable for use on all types of hard surfaces in cleanrooms and controlled environments, isolators, LAFs equipment exteriors and production vessels. The biocide blend is 0.2 micron filtered, filled and double bagged in a Grade C (ISO Class 7) cleanroom before irradiation using a validated process at no less than 25kGy.

Ready to use product is available in 1 litre trigger sprays and 5 litre capped bottles. For larger areas 5 litre bottles of concentrate, for dilution with water of suitable quality, are available.

Activity

Broad spectrum activity including:

Bactericidal Activity	<i>P.aeruginosa, E.coli, E.hirae, S.aureus</i>
Fungicidal Activity	<i>C.albicans</i>

Protocol

Concentrate - 5 litres of concentrated product should be diluted as a 1:3 dilution with water of suitable quality to make 20 litres of diluted disinfectant. Mix well before use.

Use biocides safely. Always read the label and product information before use.

Validation

Manufactured by an EN ISO 9001:2008, EN ISO 13485:2012 and European Medical Device Directive 93/42/EEC accredited company.

Passes EN test 1276 for bactericidal efficacy and EN 1650 for fungicidal efficacy.

Certificates of analysis and irradiation with every batch.

2 year shelf life.

Full supporting documentation is available on this product.

Product Codes

CODE	OLD CODE	LANGUAGES ON LABEL	DESCRIPTION	WATER QUALITY	STERILE / FILTERED	SIZE	UNIT OF SALE
3082580	3035240	GB;IT;ES	Klercide Quat / Biguanide Concentrate	DI	Sterile	5.0 litre	4 bottles x 5.0L
3079050	3035240	DE;FR;NL	Klercide Quat / Biguanide Concentrate	DI	Sterile	5.0 litre	4 bottles x 5.0L

For more information please contact info@ecolab.com or visit www.ecolabcc.com